



Installation guide: FLUICS Print for SciNote



- 1 Unbox the printer, make sure the yellow box attached with the magnets to the top of the printer. Connect the cables in the following order:

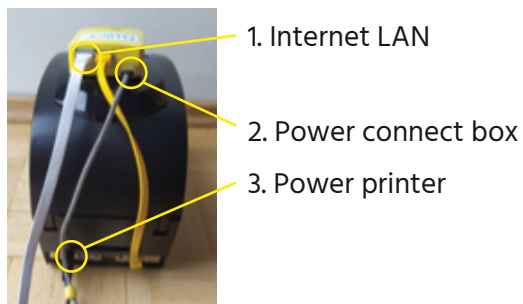

- 2 Push the "Power" button to turn on the printer and wait a few minutes. **Check that all LEDs on the yellow box are on (except for the "status" LED that blinks every 3 seconds).**

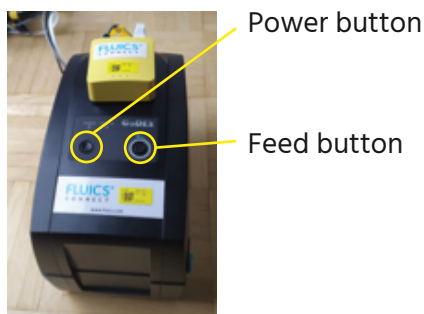

NB: Depending on your IT infrastructure it might be necessary to register the MAC address written on the bottom of the yellow box to the whitelist for devices with Internet access.

- 3 Login to your SciNote account and follow the installation steps as described on the SciNote website:

scinote.net/fluics



The API key will be provided from FLUICS via email.

Only Organization administrators can enter the API key.

- 4 Enjoy the ease of printing and tracking your QR-coded labels.

For troubleshooting your printer or ordering additional labels, please contact us at support@fluics.com.

For questions related to the SciNote print Add-on or label layouts, please contact directly SciNote support@scinote.net.

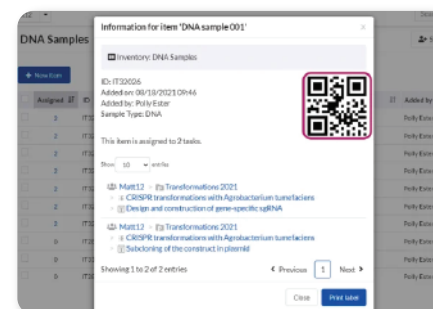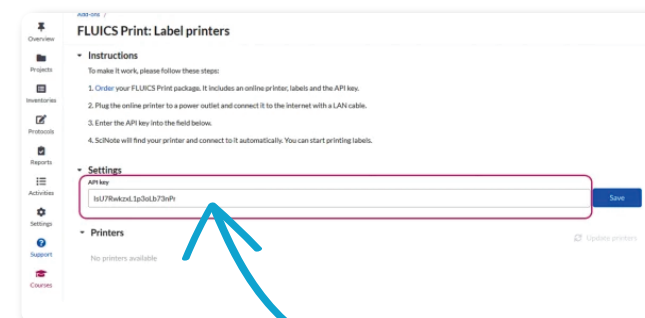

Contact details:

FLUICS GmbH
Fürstenstr. 15
80333 Munich
Germany

connect@fluics.com
+49 89 71690877

Registered in
Munich HRB 246078